

MiLab

10/07/2026 9:39 am BST

Summary

Merlin integrates with MiLab which can be accessed through the General Lab screen.

This functionality allows you to send Patients from the system onto the Lab equipment and transfer the results back to the Patient's record.

Details

Create a MiLab Request

The first step is to 'Check In' a Patient to the lab equipment.

1. Select a Patient from Reception, daylist or diary.
2. Select the 'Lab' Button.

The screenshot shows the 'Reception' screen. At the top, there's a toolbar with various icons. Below that, there's a 'Client Details' section with fields for Title, Surname, First Name, Patient, Category, Reg. Date, and Species. The 'Lab' button is highlighted with a red box.

3. Once you have entered this screen, select the 'MiLab' tab. Select the 'Create Lab Request' button to proceed.

The screenshot shows the 'General Lab' screen. At the top, there's a toolbar with various icons. Below that, there's a 'Create Lab Request' button highlighted with a red box. The 'MiLab' tab is also highlighted with a red box.

The screenshot shows the 'MiLab New Order' window. At the top, there's a toolbar with 'Create Request' and 'Exit' buttons. Below that, there's a form with fields for Surgeon, MiLab Gender, Species, and Breed. The 'Create Request' button is highlighted with a red box.

4. Once selected, the below window will display. Select the Request/s from the list and again, select the 'Create Request' button to proceed and send the request to the MiLab machine.

Results can also be viewed via the Patient's history screen.

Any lab requests done directly from the MiLab machine will still show in the results list but they will not be saved in the patients history. To do this click on the test result, select the Transfer button, search for the relevant client and patient and press transfer.

