

Abaxis Fuse

10/07/2026 9:43 am BST


Summary

The Abaxis Vet Scan Fuse integration allows the ordering of tests from within Merlin directly to the Abaxis lab equipment, when results are available they are automatically saved to the patients history.

Details


How do I create a Lab Request?

1. Select a Patient and enter the Lab Screen.

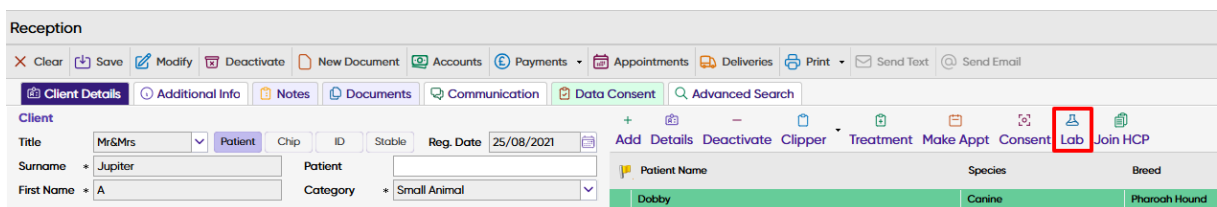


The screenshot shows the 'Reception' screen in the Merlin software. The top navigation bar includes icons for various functions. The 'Lab' icon, which represents a test tube, is highlighted with a red rectangular box. Below the navigation bar, the 'Client Details' tab is active, showing patient information such as Title, Surname, First Name, and Reg. Date.

2. Select the 'Vet Scan Fuse' tab.

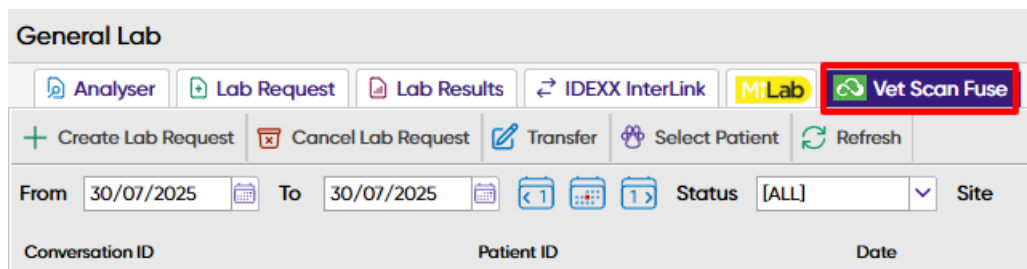

The screenshot shows the 'General Lab' screen. The 'Vet Scan Fuse' button, which features a green circular icon with a white test tube, is highlighted with a red rectangular box. Other buttons visible include 'Analyser', 'Lab Request', 'Lab Results', 'IDEXX InterLink', and 'MiLab'. Below the buttons, there are fields for 'From' and 'To' dates, a 'Status' dropdown menu, and a 'Site' dropdown menu.

3. Then, select 'Create Lab Request'.

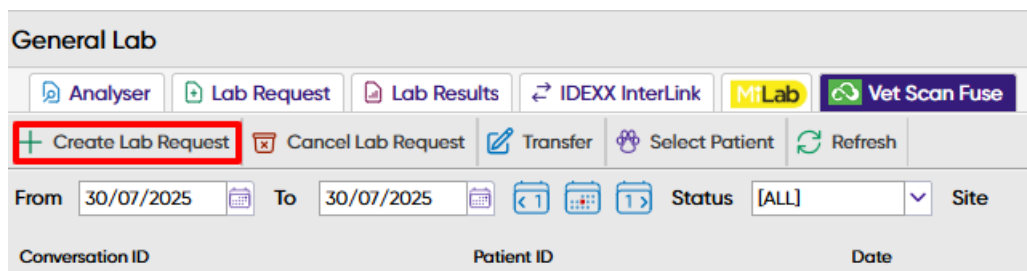

The screenshot shows the 'General Lab' screen with the 'Create Lab Request' button highlighted by a red box. This button is located in the top left of the main content area, next to the 'Analyser' button. The 'Vet Scan Fuse' button remains highlighted with a red box. The interface includes various filters and search options at the bottom.

4. The following window will appear. From this window, tick the required lab requests.

Vet Scan Fuse New Order

Create Request **Exit** Note: Hematology machines cannot accept other species

Sites:

Surgeon:

Vet Scan Gender:

Vet Scan Species:

Breed:

Lab Requests

Chemistry

☐ Avian/Reptilian Profile Plus(AR)

☐ Comprehensive Diagnostic(CDP)

☐ Critical Care Plus(CC)

☐ Equine Profile Plus(EPP)

☐ Kidney Profile Plus(KPP)

5. When selecting a request, you are prompted to select the relevant Abaxis Device.

Select the Devices

Not Specified

VetScan VS2SN: 0000V25003Status: Online

6. Once you have selected the necessary requests and selected the relevant devices, select 'Create Request' to generate.

Vet Scan Fuse New Order

Create Request

7. The lab request will display under the Vet Scan Fuse Tab. It will remain under this tab with the updated status displayed in the 'Status' column.

General Lab						
From	10/07/2026	To	10/07/2026	Status	[ALL]	Site
Conversation ID	Patient ID	Date	Client	Patient	Status	
368cd1c4-5704-4a7a-a7dd-659d045ebbf7	25502	10/07/2026 09:27			WAITING-FOR-RESPONSE	
b2c0c0a8-3fad-488f-a508-7001a3c7dec0	91253	10/07/2026 08:13			COMPLETED	

8. You can double click on the request to view details.

9. From this screen you can cancel the request or print the lab request.

Lab Results for Tony Stark

Print **Cancel Tests** **Exit**

Vet Fuse Lab Results

Patient ID: Doctor:

Patient: Conversation ID:

Results On: Status:

Species:

Result Type:


Machine Name:

Lab Requests

Chemistry

☐ Prep Profile II(Prep)

NOTE: any lab requests done directly from the VetScan machine will still show in the results list but they will not be saved in the patients history. To do this click on the test result, select the Transfer button, search for the relevant client and patient and press transfer.

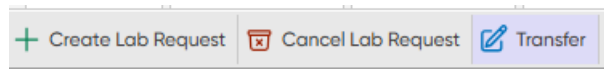

How do I view Lab Results?

Results will be displayed under the Vet Scan Fuse Tab. These can be accessed by double clicking the relevant record and selecting 'print' in the top left corner.

Alternatively, when in the patient record, select 'Patient Lab', update the date range and results can be accessed from here.

How to Setup Integration

Please contact your Territory Manager (TM) and our installations team will assist with the configuration with your equipment.
